

Specra Manager 2.0 软件安装教程

1、找到软件光盘或者安装包，如下图



2、退出杀毒软件或者清理软件等妨碍安装程序进行的软件

3、双击“Setup”进入安装程序，点击 OK，一直安装至下图。

The image shows the "InstallShield Wizard" window. The title bar says "InstallShield Wizard". The main heading is "Customer Information". Below it, it says "Please enter your information." There is a small icon of a computer with a network cable. The main area contains the following text: "Please enter your name, the name of the company for whom you work and the product serial number." Below this are three input fields: "User Name:", "Company Name:", and "Serial Number:". Each field has a red text annotation: "根据实际输入" (Enter actual input) for User Name and Company Name, and "输入光盘上的序列号或者安装包里SN文档里的序列号" (Enter the serial number on the CD or in the SN document in the installation package) for Serial Number. At the bottom, there are three buttons: "< Back", "Next >", and "Cancel".




4、选择安装位置，点“Next”，一直安装到“Finish”


5、Finish 完成后，桌面将生成“Spectra Manager”快捷方式，双击打开




6、在 Spectra Manager 窗口下，打开“Program”下的“Administrative Tools”


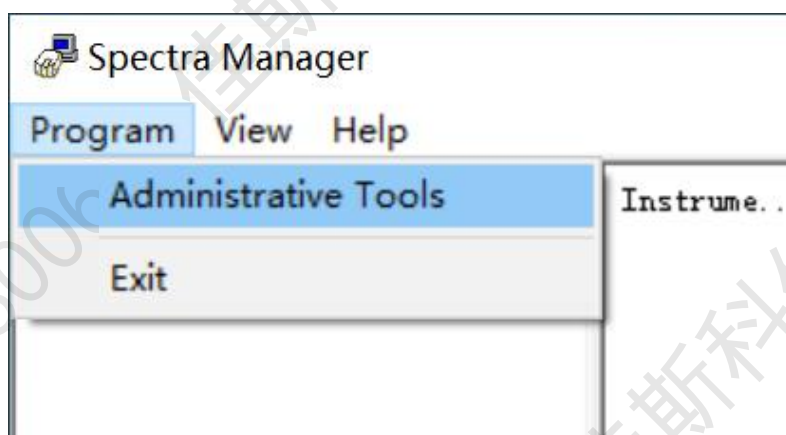


7、在 Administrative Tools 窗口下，打开“Operation”下的“Install”


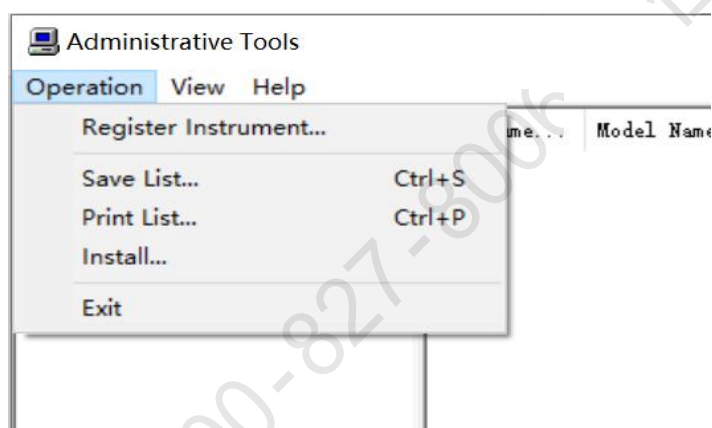


8、点击“Browse”指定驱动路径

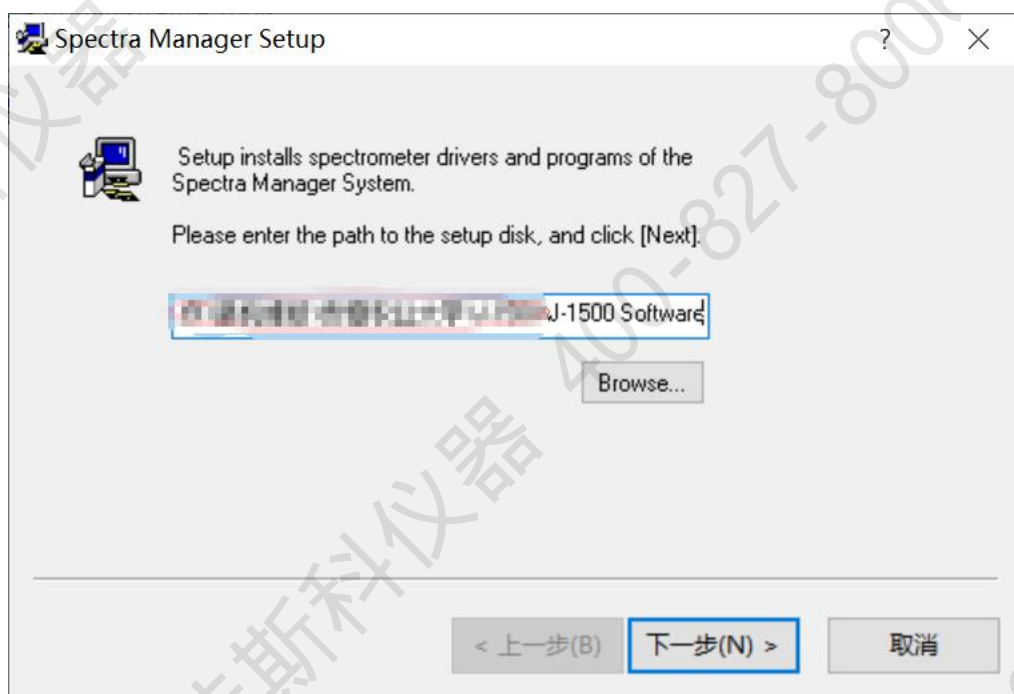

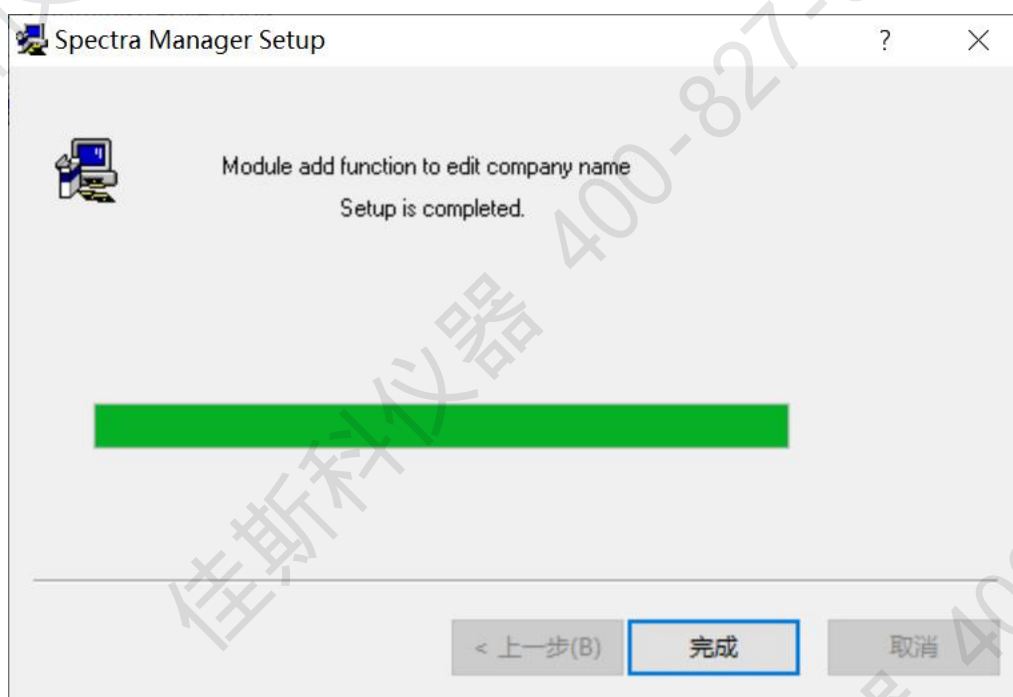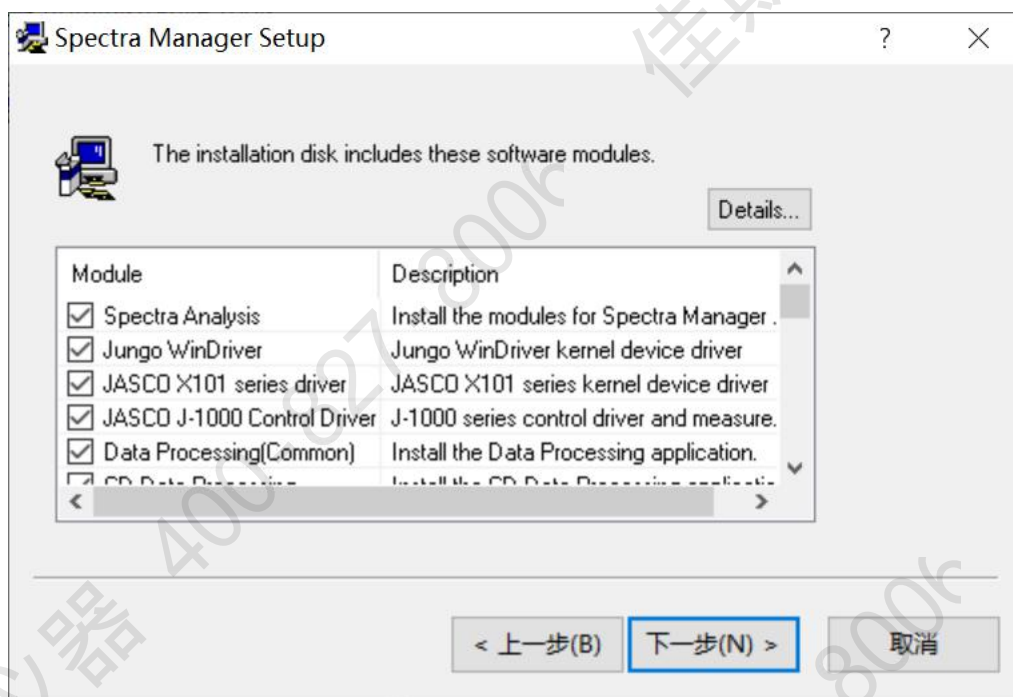

9、完成后，回到 Administrative Tools 窗口，右击 Instrument，注册仪器

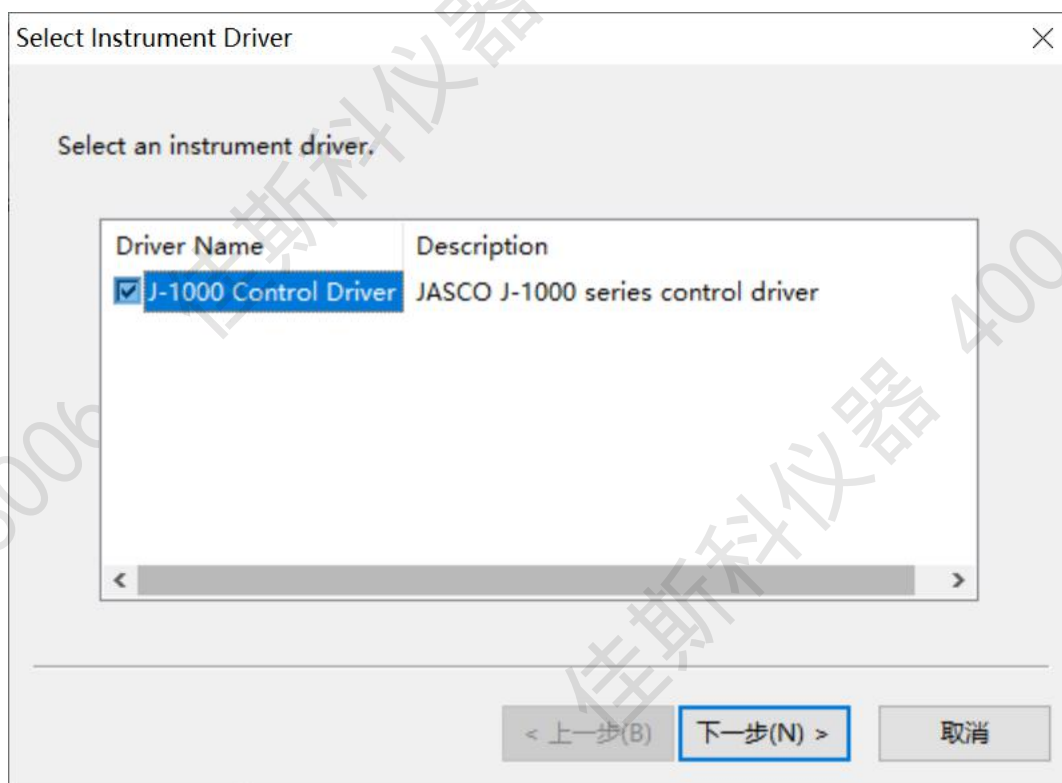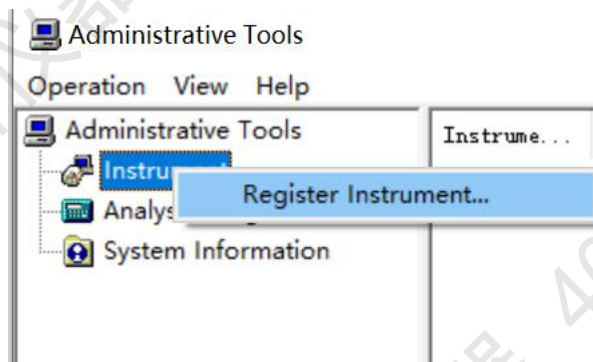

Instrument Properties


Enter an instrument name, model name, and serial number.

Instrument name:

Model name: 选择仪器的型号

Serial No.:

< 上一步(B) 完成 取消

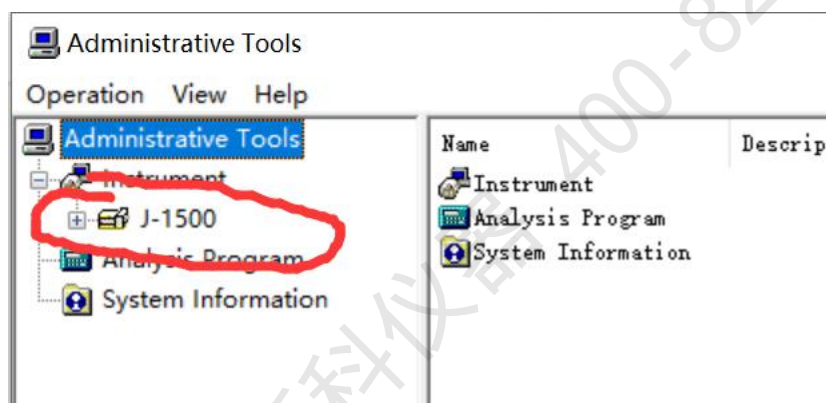

10、退出所有窗口，重新打开“Spetra Mananger”，如下图，安装完成。

